

---

# CLAB DATA-LOGGER 003

## USER'S GUIDE



CENTRE FOR MICROCOMPUTER APPLICATIONS

<http://www.cma-science.nl>

**CLAB User's Guide ver. 2.1, April 2018**

Author: Ewa Kedzierska

© CMA, Amsterdam, 2015

CMA,  
A.J. Ernststraat 169,  
1083 GT Amsterdam,  
Netherlands  
[info@cma-science.nl](mailto:info@cma-science.nl)  
[www.cma-science.nl](http://www.cma-science.nl)

CASIO and ClassPad are trademarks or registered trademarks of Casio Computer Co., Ltd.

## TABLE OF CONTENTS

<b>I. INTRODUCTION</b>	<b>5</b>
1. About CLAB	5
2. Items included with CLAB	5
3. How to use this guidebook	6
<b>II. GETTING STARTED</b>	<b>7</b>
1. CLAB overview	7
2. Power Button	8
3. Status LEDs and Speaker	9
3.1. Power LED	9
3.2. Measurement LED	10
3.3. Speaker	10
4. Powering CLAB	10
5. Memory	13
6. Sensor inputs	14
6.1. BT analog sensor plug	14
6.2. Automatic sensor detection	14
6.3. Sampling rate	15
7. Built-in Accelerometer	15
8. The ways of using CLAB	15
8.2. CLAB with a CASIO Graphic Calculator	16
8.1. CLAB with a PC computer	16
8.3. CLAB standalone	16
<b>III. USING CLAB WITH A CASIO GRAPHIC CALCULATOR AND E-CON</b>	<b>17</b>
1. CLAB and CASIO Graphic Calculators	17
2. E-CON software	17
<b>IV. USING CLAB WITH A COMPUTER AND COACH</b>	<b>18</b>
1. Coach programs	18
1.1 Coach	18
1.2 Coach Lite	18
2. Installing the CLAB driver	19
3. Typical measurement procedure in Coach	19

<b>V. USING CLAB STANDALONE</b>	<b>20</b>
3. Typical measurement procedure with CLAB standalone	20
<b>VI. INCLUDED SENSORS</b>	<b>21</b>
1. BT01 Temperature sensor	22
1. Practical information	22
2. Suggested experiments	23
3. Technical data	23
2. BT02 Voltage sensor	23
1. Practical information	24
2. Suggested experiments	24
3. Technical data	24
<b>VII. TECHNICAL SPECIFICATIONS</b>	<b>25</b>
1. Key features	25
2. Working conditions and maintenance	25
3. Software for CLAB	26
4. Warranty	26
5. Battery precautions	26
6. Safety Information	26

## I. INTRODUCTION

### 1. About CLAB

CLAB is an easy-to-use portable data-logger for Mathematics, Science and Technology Education that can be used with a CASIO Graphic Calculator equipped with the E-CON program, with a PC computer and standalone, in the field and in the classroom.

CLAB is equipped with its own processor and memory, enabling measurements at high sampling rates up to 100 000 Hz with accurate independent timing. It has FLASH memory to allow easy upgrade of the internal system software and to store measured data.

For connecting sensors CLAB has three BT (right-handed) sensor inputs, which support CMA sensors. CMA offers a wide range of sensors, which can be connected to these sensor inputs. Additionally CLAB has a built-in 3-axis accelerometer.

### 2. Items included with CLAB

The following items are included with CLAB:

- USB Power Adapter (100-240 V AC, DC 5V/1A) for powering CLAB and charging its battery
- USB cable for connecting CLAB to the computer or to the delivered USB power adapter
- BT01 Temperature sensor (-40 .. 140 °C)
- BT02 Voltage sensor (-10 .. 10 V)
- User's Quick Start Guide.



### **3. How to use this guidebook**

This guidebook is intended to provide instructions for the basic operations of the CLAB data-logger. The chapters in this guidebook include:

**II. Getting started** – provides start-up information for the learning of data-logger, and offers an overview of basic operations of CLAB.

**III. Using CLAB with a CASIO Graphic Calculator and E-CON software** – explains how to use CLAB with a CASIO Graphic Calculator.

**IV. Using CLAB with a computer and Coach software** – explains how to use CLAB with a PC computer and Coach software.

**V. Using CLAB standalone** - explains how to use CLAB standalone.


**VI. Included sensors** – provides technical specifications of included in the CLAB package sensors.

**VII. Technical specifications** – provides technical specifications, working conditions and maintenance.

This User's Guide and more on CLAB is also available on-line as a free download at CMA website [www.cma-science.nl](http://www.cma-science.nl).

## II. GETTING STARTED

### 1. CLAB overview

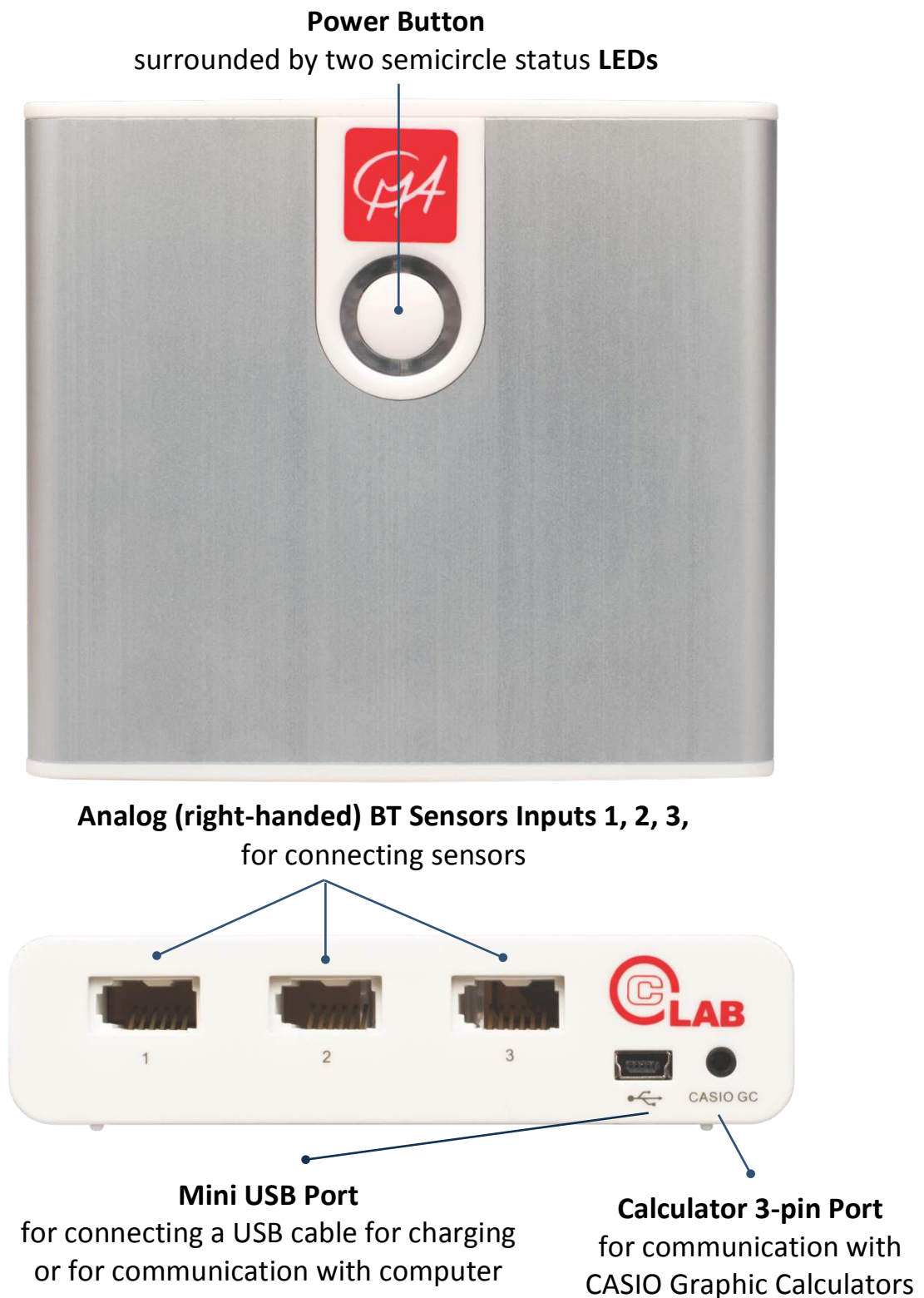

CLAB has built-in a 3-axis Accelerometer and Speaker.

## 2. Power Button

The Power Button of **CLAB** is used to:

- Turning the device on and off.
- Rebooting the device. Use this feature when **CLAB** “locks up” and does not respond to commands. Data and measurement settings may be lost.
- Handling data collection process.


The detailed operations of the Power Button:

CLAB Status	POWER BUTTON	ACTION
OFF	Press	Turns <b>CLAB</b> on The left LED blinks red and turns green when the device is ready
ON	Press and hold for 2 sec and release	Turns <b>CLAB</b> off Press until the left LED becomes red then release, the device turns off.
ON	Press and hold for 6 sec and release	Resets <b>CLAB</b> (use only when the device does not respond) Press and hold, after 2 sec the left LED becomes red, after 6 sec the LED starts to blink, then release, the rebooting starts.
ON CLAB waiting for triggering	Press	Manually triggers the data collection Press the button when the right LED is blinking blue.
ON CLAB is sampling	Press	Stops data collection Press the button when the right LED is blinking green.
ON CLAB ready for data collection	Press	Starts data collection or Collects a single measurement (Only available when the device works standalone) Press the button when the right LED is blue.



### 3. Status LEDs and Speaker

The two semicircles multicolored LEDs surrounding the Power Button together with built-in speaker provide information about the operating status of the CLAB.

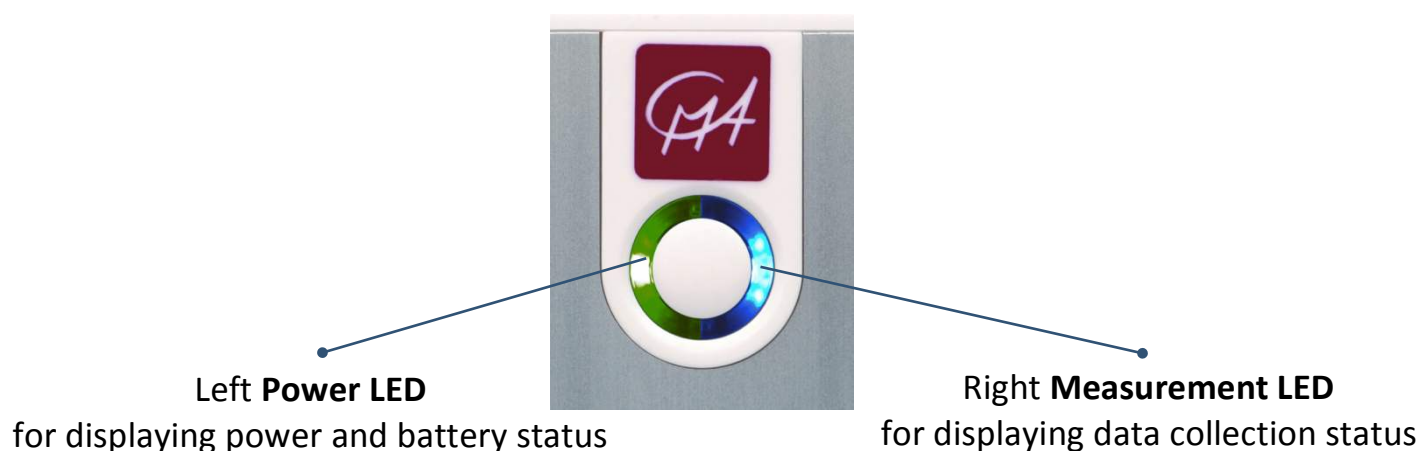

#### 3.1. Power LED

The left Power LED indicates power and battery status.

POWER LED	CLAB STATE
OFF	CLAB is turned Off and not being charged, or CLAB is in Power Save Mode
ON, together with right LED The LEDs color depend on the battery level Orange: battery is being charged Green: battery fully charged	CLAB is Off and being charged (connected to USB or to its power adapter)
ON The LED color depends on the battery level Green: battery full (battery level is between 80% and 100%) Blue: battery good (battery level is between 20% and 80%) Red: battery low (battery level is less than 20 % and needs to be charged)	CLAB is turned On and not being charged (not connected to USB or to its power adapter)  Note: powering CLAB in this situation may cause faulty battery level indication

### 3.2. Measurement LED

The right Measurement LED indicates the data-collection status.

MEASUREMENT LED	CLAB STATE
OFF	No data collection
ON Blue	Ready for data collection
ON Blinking Blue	Waiting for trigger conditions Data collection starts automatically after the trigger conditions are met. To start the data collection manually, press the Power Button.
ON Blinking Green	Sampling data
ON Green	Measurement data present in CLAB's memory
ON Red	Error

### 3.3. Speaker

Speaker generates different sounds and indicates events.

SPEAKER	CLAB STATE
Sound 1	Power On
Sound 2	Data collection starts
Sound 3	Data collection ends
Sound 4	Count down - sound n times (= number of countdown)

## 4. Powering CLAB


An internal rechargeable battery (Lithium-Ion battery, 3.7 V, 1200 mAh) located in the internal compartment, powers **CLAB**. **CLAB** is delivered with its battery partially charged. The device will work as long as the battery provides power or it is powered via the power adapter. The Power LED indicates the actual battery level when **CLAB** is turned on (see “3. Status LEDs and Speaker”).

---

#### **Note:**

- *In case **CLAB** is connected to USB or to the USB Power Adapter the Power LED will show nearly always Green.*
-

Before the battery becomes too low for operating the device the Power LED turns red to warn that the battery level is critical (lower than 20 %). If you continue working without charging, the device will shut down. A fully discharged battery requires up to 4 hours of charge time to become fully charged. But within 3 hours the battery will be charged at least up to 90 %.

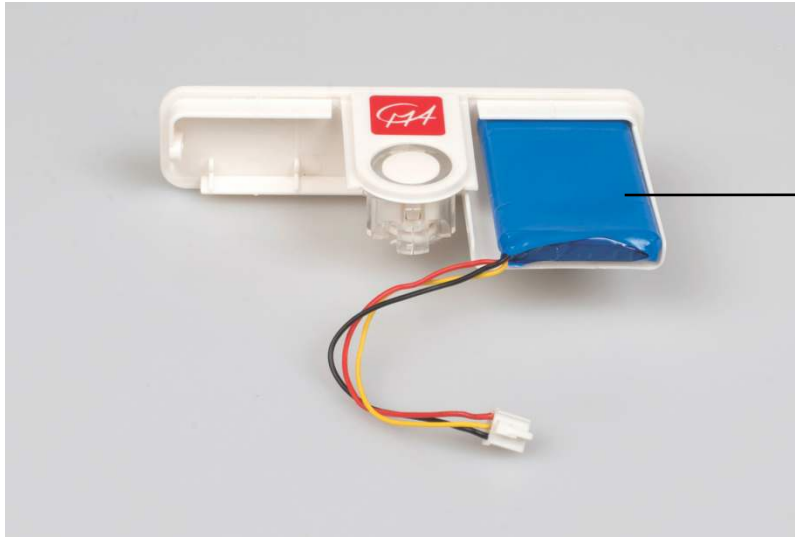

CLAB backside with rechargeable the battery placed in the battery compartment

Battery life will depend on the sensors and features used but typically you can expect to use CLAB between at least 4 hours (in case of 3 sensors connected) up to 8 hours without recharging it. To save power turn off CLAB when you do not need it. To prolong battery life, automatic power down turns off CLAB after 10 minutes of inactivity. The next time you turn it on, it is exactly as you left it.

In the case that CLAB is used standalone (see also 8.3) and the measurement settings are:

- measurement time longer than 2 hours, and
- sampling frequency less than one sample per 2 minutes,

CLAB is set into a Power-save Mode and will turn itself off between collecting data samples. It awakes 2 seconds before a new sample has to be collected. Data are saved in CLAB memory in between. In Power-save Mode the Motion detector BT55i cannot be used.

The life of the battery is more than 300 full (0 to 100%) charging cycles. If charging cycle is not full e.g. 30% to 60% it will last more charging cycles. Exposure to temperature over 35°C will significantly reduce battery life. A replacement battery can be ordered separately.

---

#### Notes:

- *CLAB can be used while the battery is being charged.*
- *CLAB typically consumes about 150 mA. If you connect many sensors then power consumption can be higher.*
- *When the USB port does not provide enough power, CLAB will not be charged or charging will be significantly slower.*

### ***To charge battery***

Use only the USB cable and USB power adapter delivered with your CLAB.

- Connect the USB cable to the mini USB port of CLAB.
- Connect the other end of the USB cable to a USB power adapter.
- Plug the USB power adapter into a standard power outlet.



- or -

- Connect the USB cable to the mini USB port of CLAB.
- Connect the other end of the USB cable into a free USB port of your computer. Notice, that the efficiency of charging via a USB port, is limited by the maximal current provided by a USB socket (500 mA for USB 2.0 and 800 mA for USB 3.0).


### ***To replace battery***

Use **only** the approved rechargeable batteries provided by CMA. To simplify the battery replacement CMA provides a complete backside of CLAB (as shown in photo) in which a new battery is already placed.

- Turn CLAB off.
- Use a small cross-slotted (Philips head) screwdriver to unscrew the screws located on the sides of CLAB.
- Very carefully move the back side of CLAB out of its housing. The battery is connected to the CLAB's PCB board via a wire and the board moves together with the plastic part.
- Very carefully disconnect the battery from the PCB by pressing the battery plug and moving it out of the battery connector on the PCB.
- Move out the Power button from the backside of CLAB. Notice that this button is made from three plastic parts, which can fall apart.
- Connect the new battery to the PCB board by inserting the battery plug into the battery connector on the board.

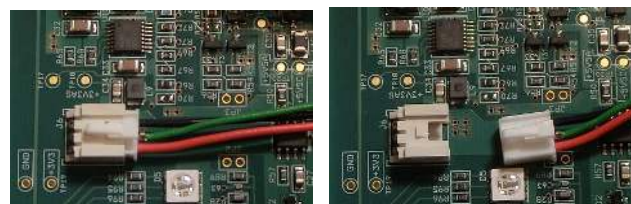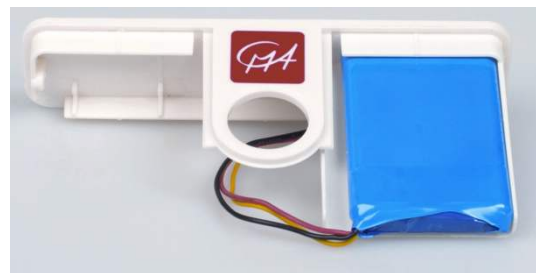

- Insert the Power button into the new backside. Take care that the parts of the Power button are placed in the correct position.
- Carefully insert the PCB board to the backside. The PCB should fit firmly.
- Keep the backside together with the board, insert the board into small rails inside the **CLAB** housing and move the board to close the housing.
- Screw the screws.
- Return your old battery for responsible recycling and do not throw it in the trash.

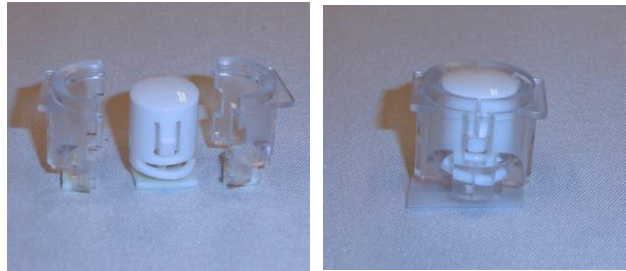

## 5. Memory

**CLAB** is equipped with 2 MB of Flash (non-volatile) memory, and 128 KB of FRAM memory.

Flash memory contains the **CLAB** internal operating system (firmware) and can be rewritten to update the firmware. As new functionality becomes available, you can download the firmware from the CMA website to your computer and update your **CLAB**. Firmware updates are also included with the Coach software.

FRAM memory of **CLAB** contains:

- the local sensor library,
- current **CLAB** experiment settings and sensor definitions,
- measured data file - which can consist of up to a total of 60,000 data points. In the Coach software it is possible to use the buffer in a cyclic way, resulting in a measurement up to 500,000 points.

If **CLAB** is switched off, all memory contents - Flash and FRAM - are maintained. FRAM is maintained as long as the battery is not completely depleted; Flash is always maintained. However, the Flash memory can be rewritten to update the local sensor library and the firmware.

### ***To update firmware***

To make use of new features for **CLAB** it is necessary to update its firmware. CMA provides a special utility program **Firmware Update**. This program can be downloaded from the CMA website or is provided with the Coach software.

The Coach program recognizes automatically when a connected to the computer **CLAB** needs to be updated and offers direct update.

## 6. Sensor inputs


For connecting sensors **CLAB** has three BT (right-handed) sensor inputs 1, 2, and 3 with input voltage ranges 0 .. 5 V and -10 .. 10 V. These inputs support all analog CMA sensors and automatically identify these sensors. Especially for **CLAB** a new CMA Ultrasonic Motion Detector BT55i, which can be connected to analog BT inputs of **CLAB**, is available.

CMA offers a wide range of sensors, for a complete list of available CMA sensors consult the CMA website. **CLAB** also supports Vernier sensors.

### 6.1. BT analog sensor plug

Sensors with BT (right-handed) plug can be connected to input channels of **CLAB**. The BT plug has 6 pins with the following connections:

Pin	Connections
1	Vin
2	Ground
3	Vres <sup>1</sup> /I2C <sup>3</sup> data
4	Auto-ID <sup>2</sup> /I2C <sup>3</sup> clock
5	5 V DC
6	Vin-low

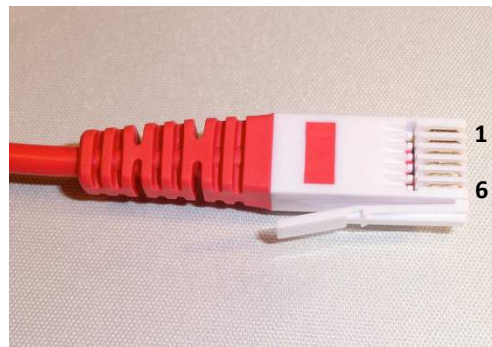

<sup>1</sup> V<sub>res</sub> - Output reference voltage (pull-up resistor 15 kΩ) for resistance measurements.

<sup>2</sup> Auto ID (pull-up resistor 10 kΩ) - Auto ID sensor detection input. The BT inputs have the possibility for automatic sensor recognition.

<sup>3</sup> For (I<sup>2</sup>C) communication between **CLAB** and intelligent sensors.

	Vin	Vin-low
Inputs	1, 2, 3	1, 2, 3
Input range	-10 .. 10 V	0 .. 5 V
Resolution (12 bit)	4.9 mV	1.2 mV
Input impedance	100 kΩ	100 kΩ

In the Coach software all **CLAB** inputs can be (simultaneously) used as counter inputs. Sensors such as the CMA Photogate or CMA Radiation sensor are by default defined as counters. Other analog sensors can be used as counters (e.g. using a Heart-rate sensor as a counter) after defining the counter conversion settings.

### 6.2. Automatic sensor detection

**CLAB** automatically detects sensors when they are connected to the sensor inputs and tries to identify the connected sensors. Most CMA sensors will be automatically identified. Some older CMA sensors or sensors connected via the

sensor adapter (e.g. CMA 4-mm to BT adapter 0519) will be not identified and need to be manually selected from the Sensor Library (on a CASIO Graphic Calculator in the E-CON software, on a PC computer in the Coach software) or a custom probe should be defined (on a CASIO Graphic Calculator in the E-CON software).

### 6.3. Sampling rate

CLAB can sample up to 3 sensors simultaneously. The sampling rate depends on the number of channels used during the data collection. For measurement via one channel maximum sampling frequency is 100 kHz, for measurement via two channels simultaneously 60 kHz and via three channels 40 kHz.

## 7. Built-in Accelerometer

CLAB has a built-in 3-axis Accelerometer, which measures acceleration in 3 ranges:  $\pm 20 \text{ m/s}^2$ ,  $\pm 40 \text{ m/s}^2$  and  $\pm 80 \text{ m/s}^2$  in the x, y and z directions. This sensor can be used in combination with other external sensors.

For the directions of x, y and z axes of the accelerometer see the image.

Holding the CLAB in the shown position should give the following values of the component accelerations:

x-component  $\approx 0 \text{ m/s}^2$

y-component  $\approx 0 \text{ m/s}^2$

z-component  $\approx -9.81 \text{ m/s}^2$ .



The exemplary activity for measurements with built-in accelerometer is delivered in Coach software (project 'Measurement with CMA CLAB' > 1.Tutorials) \*.

---

\* Starting from Coach version 6.7. Check the CMA website [www.cma-science.nl](http://www.cma-science.nl) for the latest updates.



## 8. The ways of using CLAB

CLAB can be used:

- as an interface connected to a CASIO Graphic Calculator with display of measured data on the calculator screen,
- as an interface connected to a PC computer with display of measured data on the computer screen,
- as a stand-alone data-logger.

### 8.2. CLAB with a CASIO Graphic Calculator

CLAB can be used with CASIO Graphic Calculators on which the E-CON software is installed, such as CASIO fx-CG Series, fx-9860GII Series or ClassPad II (fx-CP400). When connected to a calculator CLAB is controlled by the E-CON software running on the calculator. The collected data are transferred to the calculator and the measurement can be followed and analyzed on the calculator. To connect CLAB to a CASIO Graphic Calculator you need a data communication cable with 3-pin jack connectors (SB-62). This cable is provided with your CASIO Graphic Calculator.

### 8.1. CLAB with a PC computer

Via the provided USB cable CLAB can be connected to a USB port of a PC computer and work as a lab interface. The Coach 6 program running on the PC controls CLAB. The collected data are transferred to the computer and the measurement can be followed and analyzed on the computer screen. The USB cable is provided with your CLAB data-logger.

### 8.3. CLAB standalone

When using CLAB as a standalone device, CLAB controls the data collection process; it collects and stores the measurement data. This is very useful for remote data collection e.g. outside the school. The experimental setup is prepared using the E-CON application on a CASIO Graphic Calculator or the Coach software on a computer and uploaded to CLAB. The measured data are stored in the CLAB memory and can be downloaded into a calculator or a computer after the measurement is finished.



### III. USING CLAB WITH A CASIO GRAPHIC CALCULATOR AND E-CON

When using **CLAB** with a CASIO Graphic Calculator **CLAB** is controlled by the **E-CON** program running on the calculator. The collected data are transferred to the calculator and depending on the selected measurement mode measurement data appear during the data collection (real-time mode) or after the measurement is finished.

#### 1. CLAB and CASIO Graphic Calculators

The CASIO Graphic Calculators that have **E-CON** software installed can be used for data collection with **CLAB**.

##### *To connect CLAB to a CASIO Graphic Calculator*

Use the data communication cable with 3-pin jack connectors (SB-62) provided with your CASIO Calculator.

- Connect one end of the data communication cable to your calculator.
- Connect the other end of the cable to the CASIO Calculator port on your **CLAB**.
- Turn on **CLAB** and your CASIO Graphic Calculator.
- If you experience problems with communication make sure that both plugs are fully inserted in the communication ports.



#### 2. E-CON software

**E-CON** is a menu-driven calculator application developed by CASIO. It allows configuring experiments with **CLAB**, collecting the data via connected sensors, graphing and processing the collected data.

For more information on CASIO **E-CON** visit the following website <http://edu.casio.com/support/>.

## IV. USING CLAB WITH A COMPUTER AND COACH

When using CLAB as a lab interface connected to a PC computer CLAB stays connected to the PC via a USB port (on-line mode) and is controlled by the Coach program running on the PC. The collected data are transferred in real-time to the computer and the course of the measurement can be followed directly in the program. The recorded data are not stored in CLAB's memory and should be saved on the PC in the Coach 6 program.

### 1. Coach programs

The Coach and Coach Lite programs (version 6.6 or newer) support measurements with CLAB. Before you start to use CLAB with the computer we recommend to first install the Coach or Coach Lite program (version 6.6 or newer) on your PC.

#### 1.1 Coach<sup>1</sup>

Coach is a multimedia learning and authoring software environment for Science, Technology and Mathematics Education that offers tools to:

- collect data via interfaces and sensors
- control systems
- measure on digital videos and images
- model dynamical systems and processes
- process and analyze data.

As an authoring system Coach allows to create multimedia activities for CLAB. Exemplary Coach Activities for CLAB, prepared for use by students, are available in the Coach program.

Detailed information about the use and features of Coach 6 are described in the **Guide to Coach 6** which can be found on the CMA website ([www.cma-science.nl](http://www.cma-science.nl)).

#### 1.2 Coach Lite<sup>2</sup>

Coach Lite is a limited version of Coach, which supports measurement and control with CMA's interfaces, among others also CLAB. The data collected in measurements can be further processed; the program offers a number of data processing and data analysis tools. Exemplary Activities for CLAB, prepared for use by students, are available in Coach Lite. These Activities cannot be modified; the result of student work can be saved in Result files.

Detailed information about the use and features of Coach 6 Lite are described in the **Guide to Coach 6 Lite**, which can be found on the CMA website.

---

<sup>1</sup> Has to be purchased separate (license code needed for installation).

<sup>2</sup> Free, has to be downloaded from the CMA website [www.cma-science.nl](http://www.cma-science.nl).

## 2. Installing the CLAB driver

CLAB is equipped with HID<sup>3</sup> USB for which a standard driver is available in Windows. The driver installation is automatically executed the first time CLAB is connected to the computer<sup>4</sup>.

## 3. Typical on-line measurement procedure in Coach

*To collect data with CLAB connected to the computer<sup>5</sup>:*

- Connect CLAB to the computer via the provided USB cable.
- Turn on CLAB.
- Start Coach 6 Lite or Coach 6 and open an Activity for CLAB.
- Connect sensor(s) to the sensor input(s) on CLAB.
- CLAB automatically detects the connected sensor(s) and displays the sensor icons of the identified sensors on the CLAB screen panel. Sensors, which are not automatically identified, have to be selected manually from the Coach Sensor Library.
- The measurement in Coach is executed according to the given measurement settings. Click **Measurement Settings** to check these settings: measurement type, measuring time, sampling frequency and number of collected samples. Adjust these settings if needed.
- Follow the instructions given in the Coach Activity or directly start the measurement by clicking the green **Start** button.
- In most cases, Coach automatically starts the measurement.
  - If triggering is enabled then the measurement is started automatically when the trigger conditions are met.
  - When the type of measurement is set to **Manual** then the green **Manual Start** button appears in the Toolbar. Click this button to collect a single measurement. When keyboard input(s) is specified then the value of one or more quantities have to be typed in.
- The measurement is stopped when the specified measurement time (time-based) or specified number of samples (manual) has been reached. If you want to interrupt the measurements process, click the red **Stop** button or press **<Esc>**.
- Save your result to file via the **File** menu options **Save** or **Save As....**

---

<sup>3</sup> Human Interface Device.

<sup>4</sup> No administrator rights are needed to install this driver under all Windows versions.

<sup>5</sup> This procedure does not describe the Event-based type of measurement.

## V. USING CLAB STANDALONE

When using **CLAB** as a standalone device, **CLAB** controls the data collection process - it collects and stores the measurement data. This is very useful for remote data collection e.g. outside the school.

The experimental setup is prepared using the E-CON application on a CASIO Graphic Calculator or the Coach software on a computer and uploaded to **CLAB**. **CLAB** can be then disconnected from the calculator or computer.

The Power button of **CLAB** is used to start and stop the pre-defined experiment and LEDs and a speaker give feedback about the status of the data collection. The measured data are stored in the **CLAB** memory and can be downloaded into a calculator or a computer after the measurement is finished. Using E-CON or Coach the collected data can be further analyzed and processed.

### 1. Typical measurement procedure with **CLAB** standalone

- Prepare the experiment setting by using a CASIO Graphical Calculator with the E-CON program or a computer with the Coach program.
- Send the experiment settings to **CLAB**. The right Measurement LED of **CLAB** becomes blue when **CLAB** is ready.
- Now you can disconnect **CLAB** from your calculator or computer. If you wish (long time between preparation of the measurement and the start of the measurement) you can now turn your **CLAB** off. Settings are stored in permanent memory. After you turn **CLAB** on again the right measurement LED will show **CLAB** is ready for measurement.
- Start the measurement by pressing the Power Button.
  - If **Triggering** is enabled (the right Measurement LED is blinking blue) then the measurement is started automatically when the trigger conditions are met. To force triggering manually, press the Power Button.
  - When the type of measurement is set to **Manual** then press the Power Button to collect a single measurement.
- The right Measurement LED blinks green when **CLAB** is collecting data. If you want to stop the measurement, click the Power Button.
- For long-time measurements:
  - with time measurements longer than 2 hours, and
  - with sampling frequency of one sample per 2 minutes (or lower), and
  - without use of Motion detectors BT55i,

**CLAB** is set into a Power-save Mode and will turn itself off between collecting single data samples. It awakes 2 seconds before a new sample has to be collected. Data are saved in **CLAB** memory in between. During the measurement a short press of the Power Button will wake **CLAB** up. It will stay awake until the next sample is taken and then go back to sleep again. .

- The data collection stops after the specified measurement time (time-based mode) or specified number of samples (manual mode) has been reached.
- The Measurement LED remains green to show that measured data are stored in the **CLAB's** permanent memory. These data will stay in this memory also when you turn **CLAB** off. Only new measurement settings will cause the data to disappear.
- Turn the **CLAB** off, re-connect it to your Graphical Calculator or computer, turn it on again and retrieve the collected data into the ECON or Coach software.
- In a very rare situation it can happen that after the off-line measurement is performed **CLAB** does not start anymore. This can be caused by a problem with reading the stored data from the **CLAB's** memory. To force **CLAB** to start without reading the stored data press and hold the Power Button (3-4 sec) until the left Power LED becomes green then release the button.

## 2. Using Coach with CLAB for the standalone mode

A Coach Activity suitable for the **CLAB** standalone mode (off-line measurement)<sup>6</sup> has in its toolbar, next to the green **Start** button, two additional buttons, the **Send Settings** and **Get Data** buttons. Such activity can be used as well as for on-line measurements (**CLAB** connected to the computer during the measurement, see Chapter IV) as for off-line measurements to set up **CLAB** for remote data collection and to retrieve the collected data from its memory.

- By clicking the green **Start** button the on-line measurement is started.
- By clicking the **Send Settings** button the Experiment settings (measurement settings and sensor information) are uploaded to **CLAB**. Then **CLAB** can be disconnected from the computer and remote measurements can be performed. After the measurement is finished, **CLAB** is reconnected to the computer.
- By clicking the **Get Data** button the measurement results are downloaded to Coach. Before re-connecting **CLAB** to the computer always first turn it off. Then connect to the computer, turn it on again and open the saved Activity. Please notice that you need to be in the same Coach Activity. In the case there is no communication between Coach and **CLAB** right click the **CLAB** screen image and select **Reset Hardware**.

---

### Note:

- *Disconnecting and reconnecting **CLAB** while using it in the opened Coach Activity will give unpredictable results and can cause problems. First connect the device and then open the Activity or first (save and) close the Activity and then disconnect the device.*

---

<sup>6</sup> For off-line measurements **CLAB** does not have to be connected to the computer during the measurement. The collected data are stored in the **CLAB's** memory and have to be transferred to Coach afterwards.

## VI. INCLUDED SENSORS

Two additional sensors, Temperature and Voltage sensors are included with CLAB. Below you find technical specifications of both sensors.

### 1. BT01 Temperature sensor

The CMA BT01 Temperature sensor is a general-purpose temperature sensor that can be used to measure temperature in liquids (water, mild acidic solutions) and air. The sensor can record temperature in the range of  $-40^{\circ}\text{C}$  to  $140^{\circ}\text{C}$ . The sensing element of the sensor is a  $20\text{ k}\Omega$  NTC thermistor, which is positioned in a stainless steel tube.



The Temperature sensor is supplied calibrated in degrees Celsius  $^{\circ}\text{C}$ . The Coach 6 program allows selecting the measuring range of the sensor and shifting the sensor calibration if needed.

The sensor has a built-in resistor for automatic sensor identification.

#### 1. Practical information

The Temperature sensor can be used in a similar way as a thermometer. Here are some general guidelines for usage:

- Do not use the sensor to measure temperatures higher than  $150^{\circ}\text{C}$ , this may damage the sensor.
- Keep the sensor cable away from the heat source.
- Do not put any part of the sensor in a flame or on a hot plate.
- Avoid submerging the sensor probe beyond the stainless steel part. The handle is not waterproof.
- Always clean the sensor thoroughly after use.
- The sensor tube is constructed from stainless steel, which has a high resistance to corrosion from weak acids and alkalis. Some environments e.g. salt water may cause some discoloration of the stainless steel tube but this will not affect the sensor's performance.
- Do not use this sensor in a strong acid or strong base. A chemical reaction may cause permanent damage.
- The sensor can be left in an alkaline solution, such as NaOH, for up to 48 hours, with only minor discoloration. We do not recommend using the sensor in basic solutions which concentration is greater than 3 M.
- The maximum length of time recommended for exposure to an acid is dependent on the acid's concentration. In general we do not recommend

sensors be left to soak in acids of between 1 - 3 M concentration for longer than 48 hours. The exceptions to this are Hydrochloric acid HCL and Sulphuric acid H<sub>2</sub>SO<sub>4</sub>. The maximum exposure time for these acids are:

Acid	Maximum Exposure Time
1 M HCL	20 minutes
2 M HCL	10 minutes
3 M HCL	5 minutes
1 M H <sub>2</sub> SO <sub>4</sub>	48 hours
2 M H <sub>2</sub> SO <sub>4</sub>	20 minutes
3 M H <sub>2</sub> SO <sub>4</sub>	10 minutes

## 2. Suggested experiments

The Temperature sensor can be used in many experiments like:

- Monitoring indoor and outdoor temperatures.
- Monitoring freezing and boiling water.
- Monitoring endothermic and exothermic reactions.
- Specific heat experiments.
- Insulation investigations.
- Solar energy studies.

## 3. Technical data

<i>Temperature range</i>	-40 °C .. 140 °C
<i>Measuring element</i>	20 kΩ NTC Thermistor; non-linear
<i>Calibration (Steinhart-Hart equation)</i>	$T = [K_0 + K_1 \cdot \ln(R) + K_2 \cdot (\ln(R))^3]^{-1}$ Where T the temperature in °C, and R the measured resistance in Ω
<i>Coefficients</i>	$K_0 = 1.02119E-03$ ; $K_1 = 2.22468E-04$ ; $K_2 = 1.33342E-07$
<i>Maximum temperature that sensor can tolerate without damage</i>	150 °C
<i>Accuracy</i>	± 2°C at -40 °C ± 0.6 at 30 °C ± 1.8°C at 140 °C
<i>Response time (time for 90% change in reading)</i>	10 seconds (in water with stirring) 270 seconds (in still air) 80 seconds (in moving air)

## 2. BT02 Voltage sensor

The CMA BT02 Voltage sensor is a generic probe that can be used to measure voltage between -10 .. 10 Volts. It has two banana (4-mm) plugs for easy connection.

Note that this sensor has a direct connection to the inputs of CLAB. The sensor does not have differential inputs, and will not behave like a traditional voltmeter.



To measure a potential difference between two points in a circuit, connect the red and black plugs to the two points. A positive reading means the red plug is at a higher potential than the black plug.

The sensor has a built-in resistor for automatic sensor identification.

### 1. Practical information

The black lead is connected to the common ground of the interface. Therefore, when two or more voltage sensors are used at same time, the black leads have to be connected to the same point in the circuit.

### 2. Suggested experiments

The Voltage sensor can be used in variety experiments such as:

- Discharging of a capacitor
- Investigating battery life
- Electromagnetic induction, and many more.

### 3. Technical data

<i>Voltage range</i>	-10 .. 10 V Direct connection to CLAB via Pin 1 (-10 .. 10 V) and Pin 2 (ground) of sensor inputs.
<i>Calibration</i>	None $V_{in} (V) = V_{out} (V)$
<i>Resolution using 12 bit 5 V A/D converter</i>	4.9 mV
<i>Chemical tolerance</i>	None (air only)



## VII. TECHNICAL SPECIFICATIONS

### 1. Key features

PROCESSORS	PIC32MZ (512 kB SRAM)
MEMORY	2 MB Flash
POWER	Rechargeable battery Li-ion 3.7 V, 1200 mAh USB Power Adapter (100-240 V AC, DC 5V/1A)
STATUS INFORMATION	Two color LEDs (left Power and right Measurement status) One speaker
ADC RESOLUTION	12 bits
SAMPLING FREQUENCY	Max 100 000 Hz, via one channel Max 60 000 Hz, via two channels simultaneously Max 40 000 Hz, via three channels simultaneously
SENSOR INPUTS	Three sensor inputs, analog BT (right-handed) inputs
BUILT-IN SENSOR	3-axis Accelerometer ( $\pm 20 \text{ m/s}^2$ , $\pm 40 \text{ m/s}^2$ and $\pm 80 \text{ m/s}^2$ ), max. frequency 400 Hz
CALCULATOR CONNECTION	3-pin port for communication to CASIO Graphic Calculators
COMPUTER CONNECTION	Mini USB port also used for powering
CALCULATOR SOFTWARE	CASIO E-CON software
COMPUTER SOFTWARE	Coach 6 Lite (free) or Coach 6 (license required)
SENSORS INCLUDED	Temperature sensor and Voltage sensor

### 2. Working conditions and maintenance

For your safety and that of your equipment, follow these rules for handling, for cleaning CLAB, and for working more comfortably.

- Your **CLAB** is designed to be operated on a desk or hand-held.
- Your **CLAB** is not waterproof. Precautions must be taken to ensure that liquid cannot enter the unit.
- Do not expose **CLAB** to extreme heat or cold sources and to direct sunlight for extended periods of time.
- Operating **CLAB** outside the following ranges may affect performance:
  - Temperature: 5° to 40° C.
  - Relative humidity: 0 % to 90 % (noncondensing)
- Never force a connector into a port. When connecting a sensor, make sure that the sensor cable connector matches the sensor input, and that you have positioned the connector correctly.
- When cleaning **CLAB**, first turn it off and unplug the connected sensors and USB cable. Then clean with a soft, lint-free cloth. Avoid getting moisture in any openings. Do not spray liquid directly on **CLAB**.
- When charging the battery always use the provided USB power adapter.

- When replacing the battery, only use a **CLAB** replacement battery pack supplied by CMA. Make sure to handle it gently. Rough handling may cause the battery damage. Do not pull on wires. Pull only from the connectors at the ends of the wires. Dispose your old battery appropriately.

### 3. Software for CLAB

- **CASIO E-CON** – Data collection software installed on CASIO Graphic Calculators.
- **CMA Coach Lite<sup>\*</sup> for PC** – PC program to control **CLAB** with computer (free).
- **CMA Coach<sup>\*</sup> for PC** – PC program to control **CLAB** with computer (license needed), Author version which allows to create Activities for **CLAB** and additionally offers Data-Video, Modeling and Animations tools.

### 4. Warranty

The **CLAB** data-logger is warranted to be free from defects in materials and workmanship for a period of 24 months from the date of purchase provided that it has been used under normal laboratory conditions. This warranty does not apply if **CLAB** has been damaged by accident or misuse.

The **CLAB** battery is a consumable and is warranted to be free from defects in materials and workmanship for a period of 12 months from the date of purchase.

### 5. Battery precautions

- Do not place the battery in or near fire, on stoves or other high temperature locations such as direct sunlight. Doing so may cause the battery to generate heat, explode or ignite, or generate a toxic gas if placed in contact with fire.
- In the event the battery leaks and the fluid gets into one's eye, do not rub the eye. Rinse well with water and immediately seek medical care. If left untreated, the battery fluid could cause damage to the eye.
- Discard batteries according to local regulations.

### 6. Safety Information

**CLAB** is developed and produced in conformity with CE and FCC regulations.

Any changes made by the purchaser or user without first contacting the manufacturer will be subject to penalty under FCC regulations.

This product shall be handed over to your local community waste collection point for recycle of the products.




---

<sup>\*</sup> Version 6.6 or newer



---

*This product is to be used for educational purposes only.  
It is not appropriate for industrial, medical, research, or commercial applications.*

---